

NIST No. 2723A (Sulfur in Diesel Fuel Oil (10mg/kg))

Sheet No.: **NSX2100V-PE-016E** Petroleum chemistry
 Measurement model: **TS-2100V System** Relevant standard: **ASTM D5453**
ASC-250L/VF-210/SD-210 **JIS K 2541**

Detection method: Ultraviolet Fluorescence method

The sulfur content in diesel oil becomes sulfur oxide (SO_x) during combustion and directly affects the atmospheric air pollution as substance of concern. For this reason, it needs to be controlled at low concentration and the quantitation method complying it is required. The sulfur analysis device (**TS-2100V**) of Mitsubishi Chemical Analytech Co., Ltd. can analyze the sulfur in diesel oil quickly with accuracy.

Sample name	Sulfur in Diesel Fuel Oil (10mg/kg)																					
Analytical item	Quantitative analysis of sulfur in combustion method																					
Standard	ASTM-D5453 : standard testing method for measuring sulfur contained in carbon hydride and fuel using an ultraviolet fluorescence detector JIS K 2541 : raw petroleum and petroleum product – sulfur content testing method – ultraviolet fluorescence method																					
Analytical principle	Ultraviolet fluorescence method: Sample is burned in argon / oxygen stream and the generated sulfur dioxide is introduced to the cell of ultraviolet irradiation. The fluorescence intensity generated by ultraviolet irradiation is measured and the amount of sulfur is calculated based on the standard curve that has been created using the standard sulfur sample. $\text{Organic-S} + \text{O}_2 \rightarrow \text{SO}_2 + \text{CO}_2$ (combustion) $\text{SO}_2 + \text{h}\nu \rightarrow \text{SO}_2 + \text{h}\nu_2$ (ultraviolet fluorescence)																					
Result of sulfur analysis	<table border="1"> <thead> <tr> <th rowspan="2">Sample name</th><th colspan="5">TS-2100V analysis value (S mg/kg)</th></tr> <tr> <th>1</th><th>2</th><th>3</th><th>Average</th><th>RSD (%)</th></tr> </thead> <tbody> <tr> <td>Sulfur in Diesel Fuel Oil (10mg/kg)</td><td>10.61</td><td>10.56</td><td>10.56</td><td>10.58</td><td>0.29</td></tr> </tbody> </table>					Sample name	TS-2100V analysis value (S mg/kg)					1	2	3	Average	RSD (%)	Sulfur in Diesel Fuel Oil (10mg/kg)	10.61	10.56	10.56	10.58	0.29
Sample name	TS-2100V analysis value (S mg/kg)																					
	1	2	3	Average	RSD (%)																	
Sulfur in Diesel Fuel Oil (10mg/kg)	10.61	10.56	10.56	10.58	0.29																	
Vertical combustion method																						
Required analysis time	Pretreatment (---) minutes, Measurement (4) minutes																					
Vertical type	Total (4) minutes/ (1) measurement																					

*This sheet is provided as a reference and does not guarantee analytical values. Optimal conditions may vary depending on external factors, such as the analysis environment, and the nature of the sample.

Conditions of sulfur analysis	Measurement condition		Gas flow rate	
	Temperature of electric furnace		Ar..... 100mL/min	
Vertical combustion method	Inlet Temp	900°C	O ₂	500mL/min
	Outlet Temp	1,000°C	Ar time: 30sec	
	PMT Range	High (for Low concentration)	O ₂ time: 120sec	
	Standard sample for standard curve: S_Dibutyl disulfide / toluene			
	0, 1, 10, 50µg/mL × 30µL			
	Amount of introduced sample: 30µL			
	- The sample was not diluted, but was introduced directly.			