

Sheet No.

GT-310-ME-016E Pharmaceuticals and Cosmetics

Standardization of Potassium Permanganate Solution (Japanese Pharmacopoeia (1))

— 1/3

Reference standard: The Japanese Pharmacopoeia 18th edition

Outline

The section titled “General Tests / 9.21 Standard Solutions for Volumetric Analysis” in the Japanese Pharmacopoeia provides information on the preparation and standardization of standard solutions for volumetric analysis. The standardization of 0.02 mol/L potassium permanganate solution involves titrating sodium oxalate with potassium permanganate solution until a light red color persists for 30 seconds. In this application sheet, potentiometric titration was performed using “Return time” method to determine the end point. End point is determined correspond to the light red color persists for 30 seconds. For the common method using the inflection point as the end point, refer to Application Sheet GT310-ME-015E “Standardization of Potassium Permanganate Solution (Japanese Pharmacopoeia (2))”. The measurement results showed a relative standard deviation (RSD) of less than 1 %.

Principle

Redox titration of sodium oxalate is performed using potassium permanganate. Under acidic conditions, potassium permanganate acts as a pentavalent oxidizing agent, resulting in the reaction shown below. Since this reaction rate is slow, the solution is heated to accelerate the reaction. Since manganese (II) ions produced by the redox reaction act as a catalyst, potassium permanganate is reacted in advance to generate manganese (II) ions. Afterward, the solution is reheated, and titration is performed. Titration is performed while recording the potential difference between a platinum electrode and the reference electrode. The end point is detected when the potential corresponding to the light red color persists for 30 seconds. The factor of the 0.02 mol/L potassium permanganate solution is calculated from the mass of sodium oxalate and the volume of titrant used to reach the end point.



Apparatus

Automatic titrator: GT-310
Electrodes: PLATINUM ELECTRODE, L=105 (GTPT1B),
REFERENCE ELECTRODE, L=105 (D-J) (GTRE10B) (Outer solution: 1 mol/L potassium nitrate solution, Inner solution: 1 mol/L potassium chloride solution)

Reagents

[Titrant] ■ Potassium permanganate solution 0.02 mol/L (as specified in Japanese Pharmacopoeia)
[Reagents] ■ Sodium oxalate (certified reference material)
■ Diluted sulfuric acid (1:20): Sulfuric acid (special grade) was diluted 20 times by volume with pure water.

Sheet No.

GT-310-ME-016E

Standardization of Potassium Permanganate Solution (Japanese Pharmacopoeia (1))

2/3

Analytical Procedure

1. Sodium oxalate was dried between 150 °C to 200 °C for 1 to 1.5 hours (in this application sheet, 175 °C for 1.5 hours), then cooled in a desiccator (with silica gel).
2. Approximately 0.3 g was accurately weighed into a 500 mL beaker* and dissolved in 30 mL of purified water.
3. 250 mL of diluted sulfuric acid (1:20) was then added, and the solution was warmed to between 30 °C and 35 °C. Under gentle stirring, 40 mL of solution was quickly added (the GT-310BRT dispensing function was used). The mixture was left to stand until the red color disappeared.
4. After heating to between 55 °C and 60 °C, titration was performed using the 0.02 mol/L potassium permanganate solution.

* Although a 500 mL conical flask is used in the standard, a beaker was used instead. This was due to the small solvent volume, which would not allow full immersion of the electrode in a solution.

[Calculation]

$$\text{Factor (f)} = W / X1 / (A1 + 40) \times X2 \times 10$$

- W: Mass of sodium oxalate (g)
 X1: Mass of sodium oxalate equivalent to 1 mL of 0.02 mol/L potassium permanganate solution (= 6.7 mg/mL)
 A1: Volume of 0.02 mol/L potassium permanganate solution required for titration (mL)
 40: Volume of 0.02 mol/L potassium permanganate solution added in Step 3 (= 40 mL)
 X2: Purity of sodium oxalate (= 99.96 %)
 10: Factor to convert units, g · % to mg

Other Requirements

- Confirm reagent labels and safety data sheets for safety.
- Wear safety goggles, gloves, and/or other safety equipment when handling reagents.
- Replace the reference electrode inner and outer solutions at regular intervals.

Measurement Results


Sample	Sample amount (g)	Titration volume (mL)	Factor	Average	RSD (%)
	0.3056	5.6980	0.998		
Sodium oxalate	0.3066	5.8320	0.998	0.997	0.2
	0.3083	6.2820	0.994		

Sheet No.

GT-310-ME-016E

**Standardization of Potassium Permanganate Solution
(Japanese Pharmacopoeia (1))**

3/3

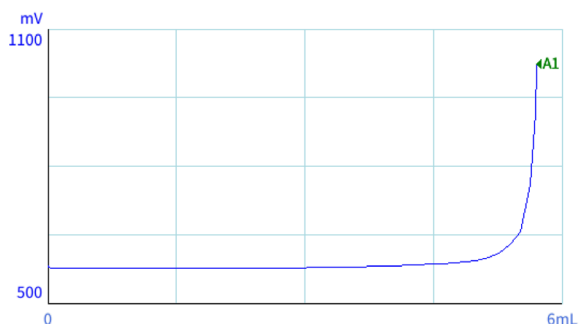

Sample name:	Sodium oxalate	
End point:	5.6980 mL	1,024.4 mV
Start of measurement:	0.000 mL	582.5 mV
End of measurement:	5.6980 mL	1024.4 mV
Measurement time:	6 min 17 s	

■ Default values were used for parameters not listed below.

Stirrer speed:	3.5
Titration mode:	Return time: R-TIME
Detector:	mV
Initial wait time:	10 s
Drop volume control:	Individual
Max. drop volume:	150 μ L
Min. drop volume:	10 μ L
Gain:	2.5
Drop cycle:	7 s
End evaluation time:	30 s
Target potential:	950 mV
Max. titration volume:	20 mL

* This application sheet is provided as reference, and does not assure the measurement results. Please consider the analysis environment, external factors and sample nature for optimal conditions before the measurement.